**English**

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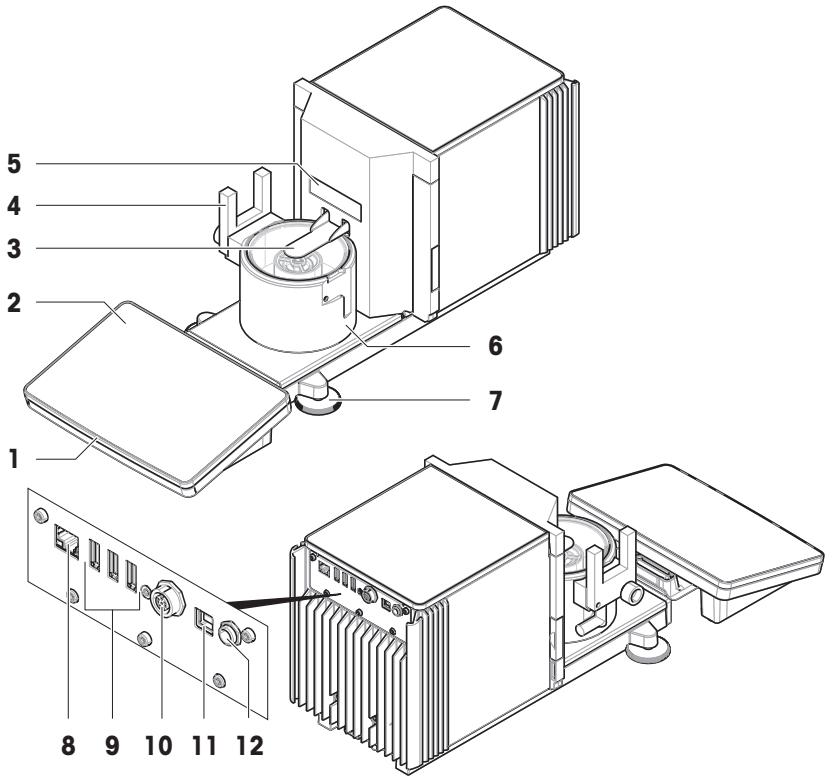
## Overview balance

en

1	StatusLight	7	Leveling feet
2	Terminal	8	Ethernet port
3	Evaporation trap door	9	USB-A ports (to device)
4	Light barrier	10	Socket for terminal connection cable
5	Balance type plate	11	USB-B port (to host)
6	Pipetting base	12	Socket for AC/DC adapter

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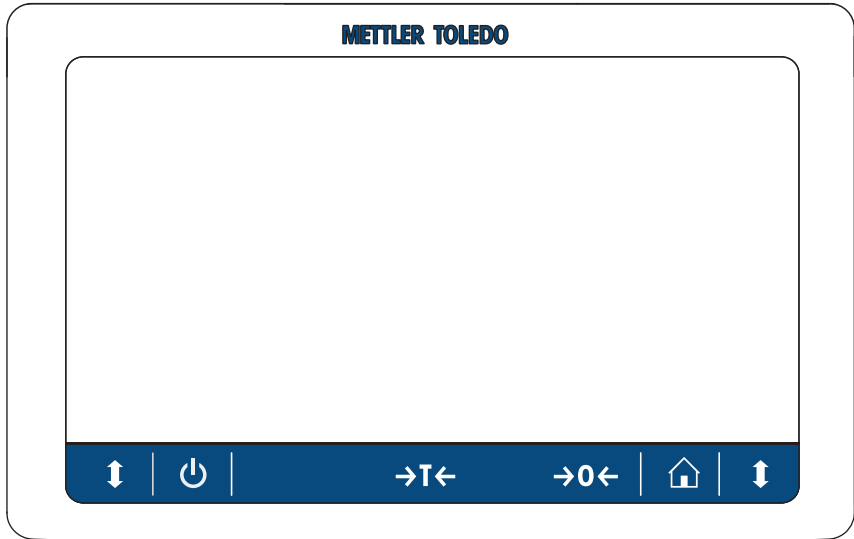
## Overview terminal

en

	Name	Description
	Standby	By tapping  , the balance is not completely switched off but goes into standby mode. To switch the balance completely off, it must be unplugged from the power supply.  <b>Note</b> Do not disconnect the balance from the power supply unless the balance is not used for an extended period of time. After switching on the instrument, it must warm up before giving accurate results.
→T←	Tare	Tares the balance.  This function is used between subsequent measurements. After taring the balance, the screen shows <b>net</b> which indicates that all displayed weights are net.
→0←	Zero	Zeroes the balance.  The balance must always be zeroed before starting the weighing process. After zeroing, the balance sets a new zero point.
	Home	To return from any menu level to the main weighing screen.
	Open/close door	Opens or closes the evaporation trap door.

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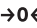
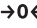
---

User Manual **Pipette Calibration Balances**

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---

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# Table of Contents

<b>1</b>	<b>Introduction</b>	<b>3</b>
1.1	Further documents and information .....	3
1.2	Acronyms and abbreviations .....	3
1.3	Compliance information .....	3
<b>2</b>	<b>Safety Information</b>	<b>4</b>
2.1	Definitions of signal words and warning symbols .....	4
2.2	Product-specific safety information .....	4
<b>3</b>	<b>Design and Function</b>	<b>5</b>
3.1	Overview .....	5
3.2	User interface .....	6
3.2.1	Main sections at a glance .....	6
3.2.2	Main weighing screen .....	6
<b>4</b>	<b>Installation and Putting into Operation</b>	<b>7</b>
4.1	Selecting the location .....	7
4.2	Unpacking the balance .....	8
4.3	Scope of delivery .....	10
4.3.1	Balance .....	10
4.3.2	Documentation .....	10
4.3.3	Calibration kit .....	10
4.4	Installation .....	11
4.4.1	Attaching the terminal .....	11
4.4.2	Assembling the balance .....	12
4.5	Putting into operation .....	13
4.5.1	Connecting the balance .....	13
4.5.2	Switching on the balance .....	14
4.5.3	Leveling the balance .....	14
4.5.4	Performing an internal adjustment .....	14
4.5.5	Performing a function check .....	14
4.5.6	Setting the balance to standby mode .....	15
4.5.7	Switching off the balance .....	15
4.6	Performing a pipetting cycle .....	15
4.6.1	Filling the evaporation trap container .....	15
4.6.2	Zeroing the balance .....	15
4.6.3	Taring the balance .....	16
4.6.4	Performing a pipetting cycle .....	16
4.6.5	Completing the weighing .....	16
4.6.6	Emptying the pipetting container .....	16
4.7	Transporting, packing and storing .....	16
4.7.1	Transporting the balance over short distances .....	16
4.7.2	Transporting the balance over long distances .....	17
4.7.3	Packing and storing .....	17
<b>5</b>	<b>Maintenance</b>	<b>17</b>
5.1	Maintenance tasks .....	17
5.2	Emptying the containers .....	18
5.2.1	Assembling the pump .....	18
5.2.2	Emptying the pipetting container .....	18
5.2.3	Emptying the evaporation trap container .....	19
5.3	Cleaning .....	20
5.3.1	Disassembling for cleaning .....	20
5.3.2	Cleaning the balance .....	20

5.3.3	Cleaning after an overflow .....	21
5.3.4	Putting into operation after cleaning.....	21
<b>6</b>	<b>Technical Data</b>	<b>21</b>
6.1	General data .....	21
<b>7</b>	<b>Disposal</b>	<b>22</b>

## 1 Introduction

Thank you for choosing a METTLER TOLEDO balance. The balance combines high performance with ease of use.

### EULA

The software in this product is licensed under the METTLER TOLEDO End User License Agreement (EULA) for Software.

► [www.mt.com/EULA](http://www.mt.com/EULA)

When using this product you agree to the terms of the EULA.

### 1.1 Further documents and information

This document is available in other languages online.

► [www.mt.com/pjpcal](http://www.mt.com/pjpcal)

Instructions for cleaning a balance: "8 Steps to a Clean Balance"

► [www.mt.com/lab-cleaning-guide](http://www.mt.com/lab-cleaning-guide)

Search for software downloads

► [www.mt.com/labweighing-software-download](http://www.mt.com/labweighing-software-download)

Search for documents

► [www.mt.com/library](http://www.mt.com/library)

For further questions, please contact your authorized METTLER TOLEDO dealer or service representative.

► [www.mt.com/contact](http://www.mt.com/contact)

### 1.2 Acronyms and abbreviations

Original term	Explanation
ASTM	American Society for Testing and Materials
EMC	Electromagnetic Compatibility
FCC	Federal Communications Commission
GWP	Good Weighing Practice
ID	Identification
LPS	Limited Power Source
MT-SICS	METTLER TOLEDO Standard Interface Command Set
NA	Not Applicable
OIML	Organisation Internationale de Métrologie Légale (International Organization of Legal Metrology)
RFID	Radio-frequency identification
RM	Reference Manual
sd	Standard deviation
SELV	Safety Extra Low Voltage
SOP	Standard Operating Procedure
SQC	Statistical Quality Control
UM	User Manual
USB	Universal Serial Bus
USP	United States Pharmacopeia

### 1.3 Compliance information

National approval documents, e.g., the FCC Supplier Declaration of Conformity, are available online and/or included in the packaging.

► <http://www.mt.com/ComplianceSearch>

For further information, consult the Reference Manual (RM).

► [www.mt.com/XPR26PC-RM](http://www.mt.com/XPR26PC-RM)

## 2 Safety Information

Two documents named "User Manual" and "Reference Manual" are available for this instrument.

- The User Manual is printed and delivered with the instrument.
- The electronic Reference Manual contains a full description of the instrument and its use.
- Keep both documents for future reference.
- Include both documents if you transfer the instrument to other parties.

Only use the instrument according to the User Manual and the Reference Manual. If you do not use the instrument according to these documents or if the instrument is modified, the safety of the instrument may be impaired and Mettler-Toledo GmbH assumes no liability.

### 2.1 Definitions of signal words and warning symbols

Safety notes contain important information on safety issues. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results. Safety notes are marked with the following signal words and warning symbols:

#### Signal words

<b>DANGER</b>	A hazardous situation with high risk, resulting in death or severe injury if not avoided.
<b>WARNING</b>	A hazardous situation with medium risk, possibly resulting in death or severe injury if not avoided.
<b>CAUTION</b>	A hazardous situation with low risk, resulting in minor or moderate injury if not avoided.
<b>NOTICE</b>	A hazardous situation with low risk, resulting in damage to the instrument, other material damage, malfunctions and erroneous results, or loss of data.

#### Warning symbols

General hazard

Notice

### 2.2 Product-specific safety information

#### Intended use

This instrument is designed to be used by trained staff. The instrument is intended for weighing purposes.

Any other type of use and operation beyond the limits of use stated by Mettler-Toledo GmbH without consent from Mettler-Toledo GmbH is considered as not intended.

#### Responsibilities of the instrument owner

The instrument owner is the person holding the legal title to the instrument and who uses the instrument or authorizes any person to use it, or the person who is deemed by law to be the operator of the instrument. The instrument owner is responsible for the safety of all users of the instrument and third parties.

Mettler-Toledo GmbH assumes that the instrument owner trains users to safely use the instrument in their workplace and deal with potential hazards. Mettler-Toledo GmbH assumes that the instrument owner provides the necessary protective gear.

## Safety notes

### **WARNING**

#### **Death or serious injury due to electric shock**

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the METTLER TOLEDO power cable and AC/DC adapter designed for your instrument.
- 2 Connect the power cable to a grounded power outlet.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and the power plug for damage and replace them if damaged.

### **NOTICE**

#### **Damage to the instrument or malfunction due to the use of unsuitable parts**

- Only use parts from METTLER TOLEDO that are intended to be used with your instrument.

A list of spare parts and accessories can be found in the Reference Manual.

## 3 Design and Function

For further information, consult the Reference Manual (RM).

► [www.mt.com/XPR26PC-RM](http://www.mt.com/XPR26PC-RM)

### 3.1 Overview

See the sections "Overview" (graphics and legend) at the very beginning of this manual.

## 3.2 User interface

### 3.2.1 Main sections at a glance

The main weighing screen (1) is the central navigation point where all the menus and settings can be found. The **Balance menu** (2), **Methods** (3) and **Protocol** (4) open when tapping the drawers along the sides of the main weighing screen.

See also

▫ Main weighing screen ▶ Page 6

### 3.2.2 Main weighing screen

	Name	Description
1	User name	Shows the name of the current user.
2	Weighing value field	Shows the current weighing value.

	Name	Description
3	Level indicator	Indicates if the balance is leveled (green) or not (red).
4	<b>Methods</b> menu	Accesses the user-defined list of methods, tests, and alignments.
5	<b>Info weight</b>	Shows the current weighing value in another unit.
6	Warning and error message area	Shows current warning and/or error messages.
7	Drawer <b>Protocol</b>	Shows the recent weighing results.
8	Sample status <b>OK</b>	Result status indicator green: indicates that the result fulfills a set of criteria. For example: <ul style="list-style-type: none"> <li>• The balance is in level.</li> <li>• The internal adjustment was performed and ok.</li> <li>• The weighing result is within the defined tolerance range (only if tolerance is defined).</li> </ul>
9	Sample status <b>Excluded</b>	Result status indicator black: indicates that the result was excluded from the protocol.
10	Sample status <b>Not OK</b>	Result status indicator red: indicates that the result criteria are not fulfilled, e.g., "The weighing result was out of the defined tolerances".
11	Button <b>Add to protocol</b>	Adds the result to the protocol. Depending on the selected method, the button can have different functions.
12	Action bar	Contains actions referring to the current task.
13	<b>Balance menu</b>	Accesses the balance properties.
14	Method information area	Contains information about the sample, method or task IDs.
15	SmartTrac	Used as a weighing aid to define a target weight with upper and lower tolerances.
16	Weighing value area	Shows the results of the current weighing process.
17	<b>Method name</b>	Shows the name of the current method.

## 4 Installation and Putting into Operation

### 4.1 Selecting the location

A balance is a sensitive precision instrument. The location where it is placed will have a profound effect on the accuracy of the weighing results.

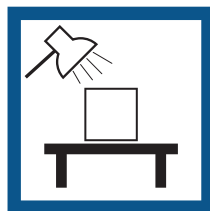
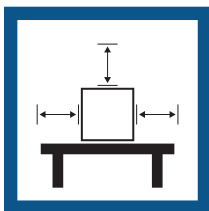
#### Requirements of the location

Place indoors on stable table

Ensure sufficient spacing

Level the instrument

Provide adequate lighting

Avoid direct sunlight

Avoid vibrations

Avoid strong drafts

Avoid temperature fluctuations

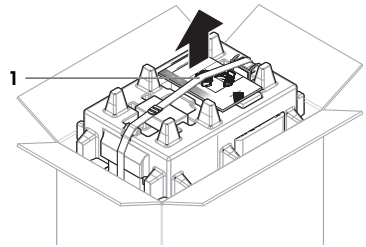
Sufficient spacing for balances: > 15 cm all around the instrument

Take into account the environmental conditions. See "Technical Data".

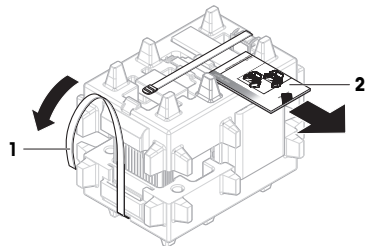
## 4.2 Unpacking the balance

Check the package, the packaging elements and the delivered components for damages. If any components are damaged, please contact your METTLER TOLEDO service representative.

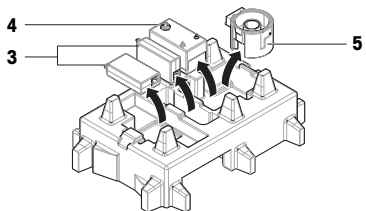
- 1 Open the box and lift the package out using the lifting strap (1).

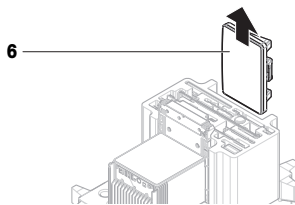
- 2 Open the lifting strap (1) and remove the User Manual (2).

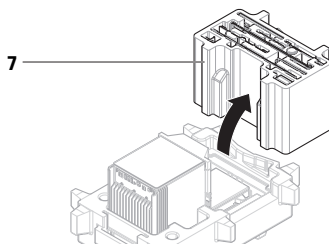
- 3 Remove the upper part of the package and remove the AC/DC adapters (3) and power cables, the pump (4), and the pipetting base with the evaporation trap container (5).

4 Carefully remove the terminal (6).

5 Carefully remove the package set with the calibration kit and other small items (7).

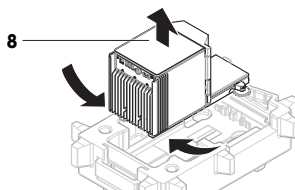
6 Carefully remove the weighing unit (8) from the bottom packaging.

7 Remove the protective bag.

8 Store all parts of the packaging in a safe place for future use.

➔ The weighing unit is ready for assembling.

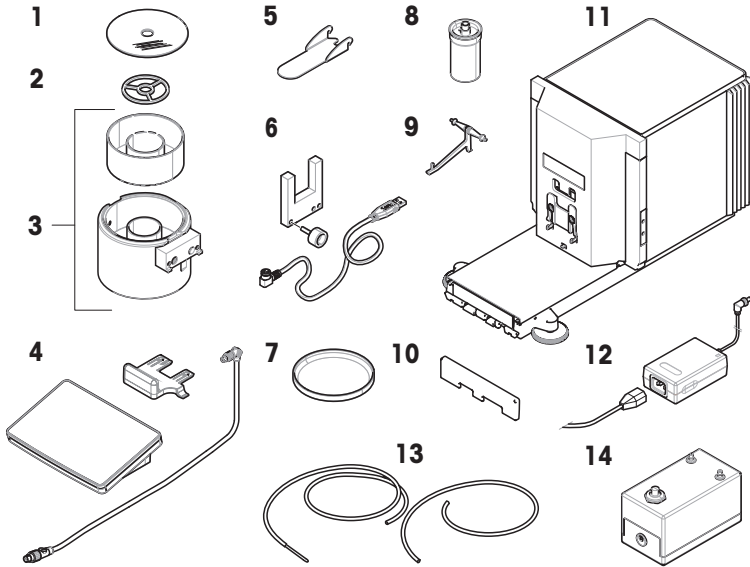
**See also**

📄 Transporting, packing and storing ▶ Page 16

### 4.3 Scope of delivery

#### 4.3.1 Balance

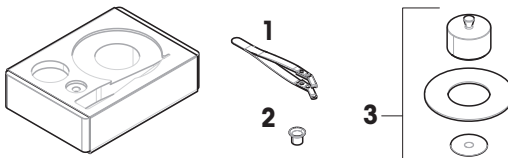
1	Evaporation trap cover	7	Drip tray
2	Centering ring	8	Pipetting container and tube
3	Pipetting base with light barrier spacer and evaporation trap container	9	Support for pipetting container
4	Terminal with terminal holder and connection cable	10	Front cover
5	Evaporation trap door	11	Weighing unit
6	Light barrier, fixing screw and cable	12	AC/DC adapter with country-specific power cable, 2 pcs
13	Pump tubes, 2 pcs (one tube with probe)	14	Pump

#### 4.3.2 Documentation

- User Manual
- Declaration of Conformity
- Production certificate

#### 4.3.3 Calibration kit

1	Tweezers	3	Weighing pan, large, with cover ring and draft shield
2	Weighing pan, small		

## 4.4 Installation

### 4.4.1 Attaching the terminal

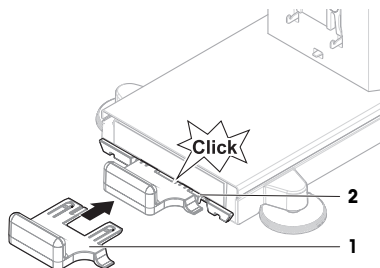
#### NOTICE

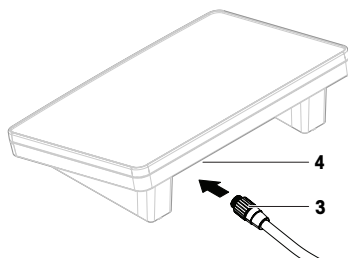
##### Damage to the cables due to careless handling

- Do not kink or twist the cables.

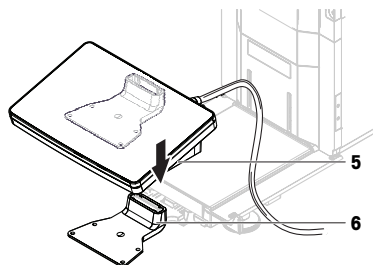
- 1 Insert the slides of the display holder (1) into the front of the weighing unit (2).

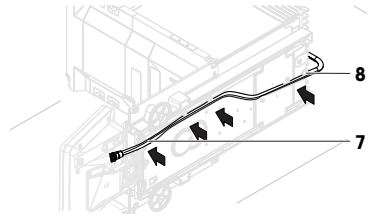
- 2 Connect the terminal cable (3) with the terminal (4). Consider the pin assignment.

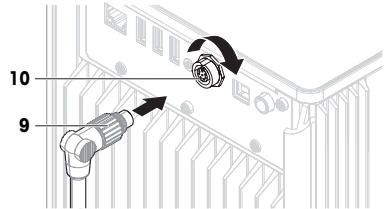
- 3 Place the terminal (5) onto the terminal holder (6).

- 4 Carefully tilt the balance to its side.
- 5 Lead the cable (7) through the cable channel (8).
- 6 Carefully put the balance back on its feet.

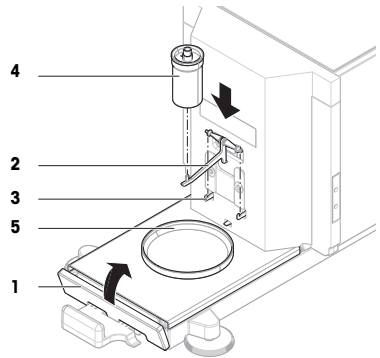
- 7 Insert the terminal cable (9) into the socket of the balance (10). Consider the pin assignment.
- ➔ The terminal is ready.

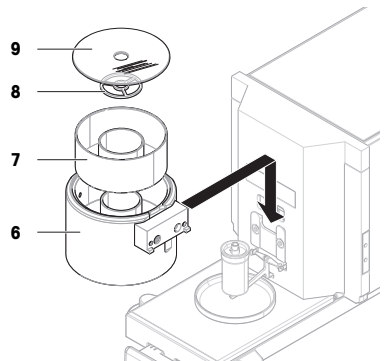
#### 4.4.2 Assembling the balance

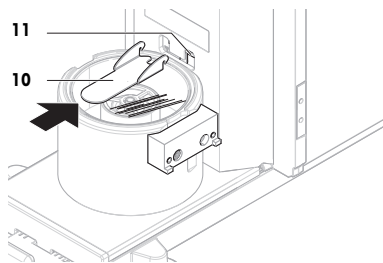
- 1 Install the front cover (1).
- 2 Carefully hang the pipetting container support (2) on the hooks (3).
  - ➔ The pipetting container support (2) hangs freely.
- 3 Carefully place the pipetting container (4) onto the pipetting container support (2).
- 4 Center the drip tray (5) underneath the pipetting container (4).

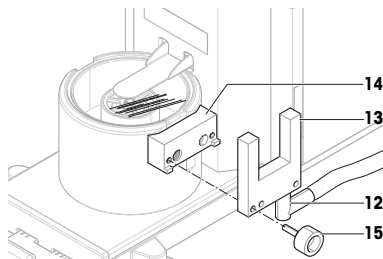
- 5 Carefully slide the pipetting base (6) with the evaporation trap container (7) vertically along the guiding rails and place it onto the drip tray.
  - ➔ The pipetting base centers the drip tray. It does not touch the pipetting container (5) or the pipetting container support (3).
- 6 Place the centering ring (8) onto the evaporation trap container (7).
- 7 Carefully place the evaporation trap cover (9) onto the evaporation trap container (7).

- 8 Slide the evaporation trap door (10) towards the back until it snaps in on the rod (11).

- 9 Connect the light barrier cable (12) to the light barrier (13).
- 10 Install the light barrier (13) on the light barrier spacer (14) using the fixing screw (15).

**Note**

By default, the light barrier and the light barrier spacer are installed on the right side of the pipetting base. Alternatively, these items can be installed on the left side of the pipetting base.

- 11 Guide the light barrier cable (12) along the side of the balance, below the side cover.
- 12 Connect the light barrier cable (12) to one of the USB-A ports on the back side of the balance.
- ➔ The balance is assembled and ready to be put into operation.

## 4.5 Putting into operation

### 4.5.1 Connecting the balance

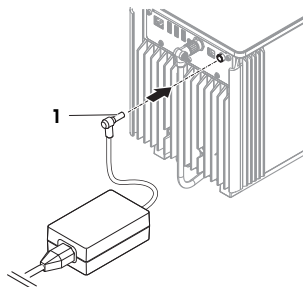
#### **WARNING**

##### **Death or serious injury due to electric shock**

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the METTLER TOLEDO power cable and AC/DC adapter designed for your instrument.
- 2 Connect the power cable to a grounded power outlet.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and the power plug for damage and replace them if damaged.

- 1 Install the cables in such a way that they cannot be damaged or interfere with operation.
- 2 Insert the plug of the AC/DC adapter (1) into the power socket of the instrument.
- 3 Secure the plug by firmly tightening the knurled nut.
- 4 Insert the plug of the power cable into a grounded power outlet that is easily accessible.

**Note**

Do not connect the instrument to a power outlet controlled by a switch. After switching on the instrument, it must warm up before giving accurate results.

## See also

📄 General data ▶ Page 21

### 4.5.2 Switching on the balance

#### EULA (End User License Agreement)

When the balance is switched on the first time, the EULA (End User License Agreement) appears on the screen.

- 1 Read the conditions.
- 2 Tap **I accept the terms in the license agreement.** and confirm with **✓ OK.**

#### Warming up

Before the balance gives reliable results, it must warm up. This takes at least 120 minutes after connecting the balance. When the balance is switched on from standby, it is ready immediately.

- The balance has warmed up.
- Press .
  - ➔ The main weighing screen appears.

When the balance is switched on, the main weighing screen appears. The display will always show the screen of the method last used before switching it off.

### 4.5.3 Leveling the balance

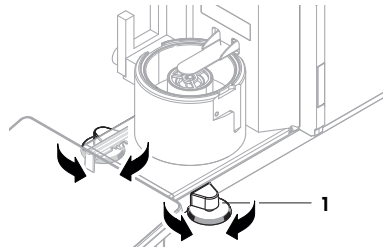
Exact horizontal and stable positioning are essential for repeatable and accurate weighing results.

If the message **Balance is out of level** appears:

- 1 Tap **▶ Level the balance.**
  - ➔ The **Leveling aid** opens.
- 2 Turn both leveling feet (**1**) as instructed on the display until the dot is in the center of the level indicator.

The leveling aid can also be accessed through the **Balance menu**:

**Navigation:** ▶ **Balance menu** >  **Leveling aid**

### 4.5.4 Performing an internal adjustment

- The adjustment **Strategy** is set to **Internal adjustment.**
- 1 Open the **Methods** section, tap , select the adjustment, and tap **▶ Start**
    - or -
    - from the main weighing screen, tap **⋮ More** and tap **Start adjustment.**
      - ➔ **Internal adjustment** is being executed.
      - ➔ When the adjustment has been completed, an overview of the adjustment results appears.
  - 2 Tap  **Print** if you want to print the results.
  - 3 Tap **✓ Finish adjustment.**
    - ➔ The balance is ready.

### 4.5.5 Performing a function check

To perform a function check, use the calibration kit delivered with the balance.

For further information, consult the Reference Manual (RM).

▶ [www.mt.com/XPR26PC-RM](http://www.mt.com/XPR26PC-RM)

#### 4.5.6 Setting the balance to standby mode

The balance can be set to standby mode by holding **⏻**. Exit the standby mode by holding **⏻** again.

#### 4.5.7 Switching off the balance

To completely switch off the balance, it must be disconnected from the power supply. By holding **⏻**, the balance goes only into standby mode.

#### Note

When the balance was completely switched off for some time, it must warm up before it can be used.

#### See also

Switching on the balance ▶ Page 14

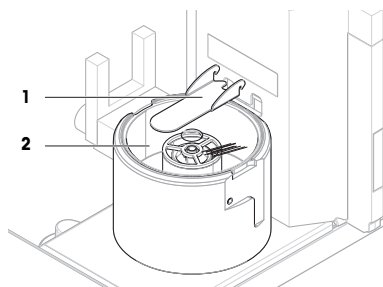
### 4.6 Performing a pipetting cycle

To calibrate the pipette, distilled water must be used. The volume of the liquid is checked by gravimetric testing.

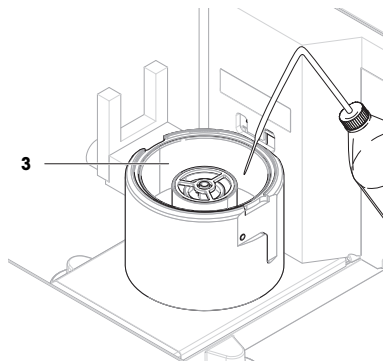
#### 4.6.1 Filling the evaporation trap container

The evaporation trap includes a container that should be filled with water. When the evaporation trap is closed, the humidity inside increases until the air is saturated. This prevents evaporation of the liquid in the pipetting container and ensures accurate weighing results.

- 1 Remove the evaporation trap door (1).
- 2 Carefully remove the evaporation trap cover (2).

- 3 Fill the evaporation trap container (3) with distilled water.
  - ➔ The evaporation trap container is at least half full (maximum 70 ml).
- 4 Reassemble the balance.
  - ➔ The evaporation trap door (1) is closed.
- 5 Wait at least **120 minutes** to allow the air in the evaporation trap to reach saturation.

#### 4.6.2 Zeroing the balance

- Press **→0←** to zero the balance.
- ➔ The balance is zeroed.

### 4.6.3 Taring the balance

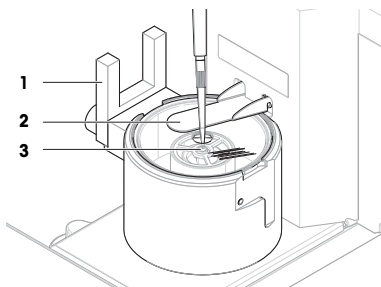
Before a subsequent measurement, the balance must be tared.

- Press →**T**← to tare the balance.
- ➔ The balance is tared. The icon **Net** appears.

### 4.6.4 Performing a pipetting cycle

- The pipette to be calibrated is prepared and set to the target volume.

- 1 Use the pipette to aspirate the target volume of distilled water.
- 2 Activate the light barrier (1), or press ↓.
  - ➔ The evaporation trap door (2) opens.
- 3 Dispense the liquid into the pipetting container (3).
- 4 Activate the light barrier (1), or press ↓.
  - ➔ The evaporation trap door (2) closes.
  - ➔ The weighing result is displayed.
- 5 Tap **+** **Add to protocol** if you want to report the weighing result.
  - ➔ The weight value is listed in the **Protocol**.

### 4.6.5 Completing the weighing

- 1 To save the **Protocol**, tap **Complete**.
  - ➔ The window **Complete task** opens.
- 2 Select an option to save or print the **Protocol**.
  - ➔ The respective dialog opens.
- 3 Follow the instructions from the wizard.
- 4 Tap **✓ Complete**.
  - ➔ The **Protocol** is saved/printed and then cleared.

### 4.6.6 Emptying the pipetting container

Regularly empty the pipetting container to prevent an overflow.

#### See also

- 📖 Emptying the pipetting container ▶ Page 18
- 📖 Cleaning after an overflow ▶ Page 21

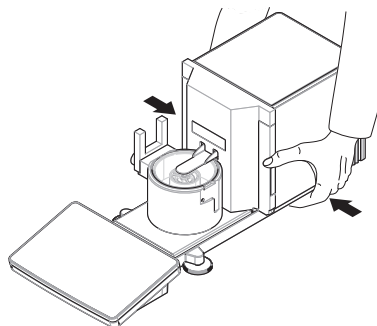
## 4.7 Transporting, packing and storing

### 4.7.1 Transporting the balance over short distances

- 1 Disconnect the AC/DC adapter and unplug all interface cables.
- 2 Hold the weighing platform with both hands and carry the balance in horizontal position to the target location.
  - Consider the requirements of the location.

If you want put the balance into operation, proceed as follows:

- 1 Connect in reverse order.
- 2 Give the balance sufficient time to warm up.
- 3 Level the balance.
- 4 Perform an internal adjustment.

**See also**

-  Selecting the location ▶ Page 7
-  Switching on the balance ▶ Page 14
-  Leveling the balance ▶ Page 14
-  Performing an internal adjustment ▶ Page 14

**4.7.2 Transporting the balance over long distances**

METTLER TOLEDO recommends using the original packaging for transportation or shipment of the balance or balance components over long distances. The elements of the original packaging are developed specifically for the balance and its components and ensure maximum protection during transportation.

**See also**

-  Unpacking the balance ▶ Page 8

**4.7.3 Packing and storing****Packing the balance**

Store all parts of packaging in a safe place. The elements of the original packaging are developed specifically for the balance and its components, and ensures maximum protection during transportation and storage.

**Storing the balance**

Only store the balance under the following conditions:

- Indoor and in the original packaging
- According to the environmental conditions, see "Technical Data"

**Note**

When storing for longer than 6 months, the rechargeable battery may become empty (only date and time get lost).

**See also**

-  Technical Data ▶ Page 21

**5 Maintenance**

To guarantee the functionality of the balance and the accuracy of the weighing results, a number of maintenance actions must be performed by the user.

For further information, consult the Reference Manual (RM).

▶ [www.mt.com/XPR26PC-RM](http://www.mt.com/XPR26PC-RM)

**5.1 Maintenance tasks**

Maintenance action	Recommended interval	Remarks
Performing an internal adjustment	<ul style="list-style-type: none"> <li>• Daily</li> <li>• After cleaning</li> <li>• After leveling</li> <li>• After changing the location</li> </ul>	see "Performing an internal adjustment"
Performing routine tests (eccentricity test, repeatability test, sensitivity test). METTLER TOLEDO recommends to at least perform a sensitivity test.	<ul style="list-style-type: none"> <li>• After cleaning</li> <li>• After assembling the balance</li> <li>• After a software update</li> <li>• Depending on your internal regulations (SOP)</li> </ul>	see "Tests" in the Reference Manual

Maintenance action	Recommended interval	Remarks
Emptying the containers	<ul style="list-style-type: none"> <li>Depending on the frequency of usage of the instrument</li> <li>Depending on your internal regulations (SOP)</li> </ul>	see "Emptying the containers"
Cleaning	<ul style="list-style-type: none"> <li>After every use</li> <li>Depending on the degree of pollution</li> <li>Depending on your internal regulations (SOP)</li> </ul>	see "Cleaning"
Updating the software	<ul style="list-style-type: none"> <li>Depending on your internal regulations (SOP).</li> <li>After a new software release.</li> </ul>	see "Software update" in the Reference Manual

### See also

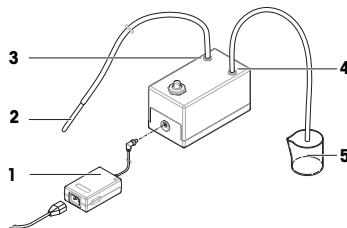
- 📄 Performing an internal adjustment ▶ Page 14
- 📄 Cleaning ▶ Page 20
- 📄 Emptying the containers ▶ Page 18

## 5.2 Emptying the containers

### 5.2.1 Assembling the pump

During normal usage, the pipetting container and the evaporation trap container are filled with water. The pump serves to remove the water in a safe and efficient manner, without having to disassemble the instrument. The inlet tube has a probe attached for easier handling. The outlet tube serves to dispose of the aspirated liquid.

- 1 Install the cables in such a way that they cannot be damaged or interfere with operation.
- 2 Insert the plug of the AC/DC adapter (1) into the power socket.
- 3 Secure the plug by firmly tightening the knurled nut.
- 4 Insert the plug of the power cable into a grounded power outlet that is easily accessible.
- 5 Connect the pump tube with the probe (2) to the nozzle IN (3).
- 6 Connect the other pump tube to the nozzle OUT (4).
- 7 Place the free end of the tube into a container suitable for liquid waste (5).

### 5.2.2 Emptying the pipetting container

The pipetting container must be regularly emptied to prevent an overflow.

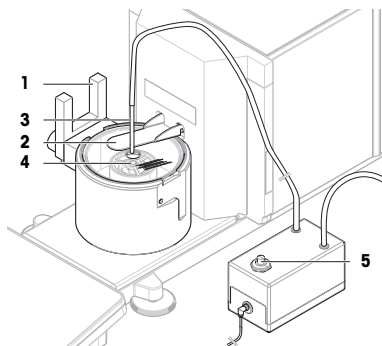
#### NOTICE

##### Damage to the pump due to inappropriate handling

- If the pump is switched on when it is not in use, it can get damaged.
- Switch off the pump when it is not in use.

- The pump is assembled.
  - The outlet pump tube is placed in a container suitable for liquid waste.
- 1 Activate the light barrier (1), or press .
    - ➔ The evaporation trap door (2) opens.
  - 2 Introduce the pump probe (3) into the pipetting container (4).
  - 3 Switch on the pump (5).
    - ➔ The liquid is pumped out and collected in the container for liquid waste.
  - 4 Switch off the pump (5).
  - 5 Reassemble the balance.
    - ➔ The balance is ready to be used.

#### See also

-  Assembling the pump ▶ Page 18

### 5.2.3 Emptying the evaporation trap container

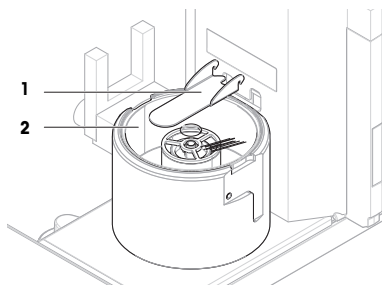
#### NOTICE

##### Damage to the pump due to inappropriate handling

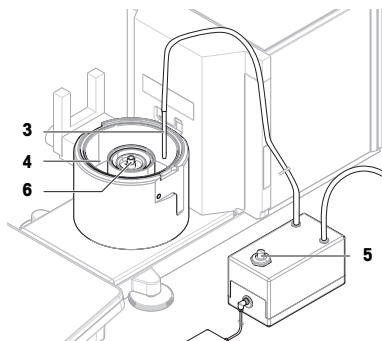
If the pump is switched on when it is not in use, it can get damaged.

- Switch off the pump when it is not in use.

- The pump is assembled.
  - The outlet pump tube is placed in a container suitable for liquid waste.
- 1 Remove the evaporation trap door (1).
  - 2 Carefully remove the evaporation trap cover (2).

- 3 Introduce the pump probe (3) into the evaporation trap container (4).
- 4 Switch on the pump (5).
  - ➔ The liquid is pumped out and collected in the container for liquid waste.
- 5 If required, use the pump to empty the pipetting container (6), see [Emptying the pipetting container ▶ Page 18].
- 6 Switch off the pump (5).
- 7 Alternatively, remove the evaporation trap container (4) to empty it.
- 8 Reassemble the balance.

#### See also

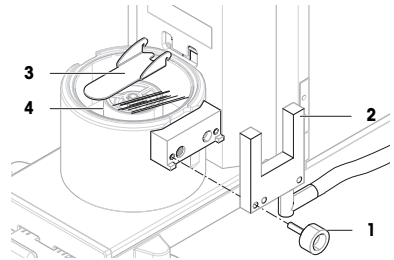
-  Assembling the pump ▶ Page 18

## 5.3 Cleaning

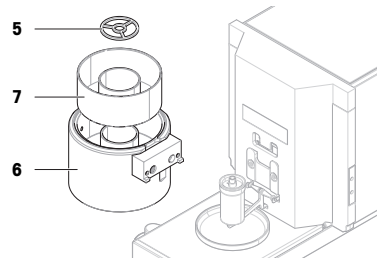
### 5.3.1 Disassembling for cleaning

- The containers are empty, see [Emptying the containers ▶ Page 18].
- The balance is switched off, see [Switching off the balance ▶ Page 15].

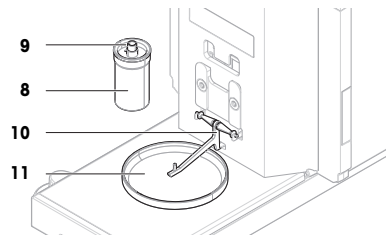
- 1 Remove the fixing screw (1) and put the light barrier (2) aside.
- 2 Remove the evaporation trap door (3).
- 3 Carefully remove the evaporation trap cover (4).

- 4 Remove the centering ring (5).
- 5 Carefully lift and remove the pipetting base (6) with the evaporation trap container (7).

- 6 Remove the pipetting container (8) and unscrew the cover holding the pipetting tube (9).
- 7 Carefully remove the pipetting container support (10).
- 8 Remove the drip tray (11).

### 5.3.2 Cleaning the balance

#### NOTICE

##### Damage to the instrument due to inappropriate cleaning methods

If liquid enters the housing, it can damage the instrument. The surface of the instrument can be damaged by certain cleaning agents, solvents, or abrasives.

- 1 Do not spray or pour liquid on the instrument.
- 2 Only use the cleaning agents specified in the Reference Manual (RM) of the instrument or the guide "8 Steps to a Clean Balance".
- 3 Only use a lightly moistened, lint-free cloth or a tissue to clean the instrument.
- 4 Wipe off any spills immediately.

For further information on cleaning a balance, consult "8 Steps to a Clean Balance".

**Cleaning around the balance**

- Remove any dirt or dust around the balance and avoid further contaminations.

**Cleaning the terminal**

- Clean the terminal with a damp cloth or a tissue and a mild cleaning agent.

**Cleaning the removable parts**

- Clean the removed part with a damp cloth or a tissue and a mild cleaning agent.

**Cleaning the weighing unit**

- 1 Disconnect the balance from the AC/DC adapter.
- 2 Use a lint-free cloth moistened with a mild cleaning agent to clean the surface of the balance.
- 3 Remove powder or dust with a disposable tissue first.
- 4 Remove sticky substances with a damp lint-free cloth and a mild solvent, e.g., isopropanol or ethanol 70%.

**5.3.3 Cleaning after an overflow**

In the event of an overflow, for example when overfilling the pipetting container, excess liquid must be removed immediately.

- 1 Use the pump to empty the evaporation trap container, see [Emptying the evaporation trap container ▶ Page 19].
- 2 Use the pump to empty the pipetting container, see [Emptying the pipetting container ▶ Page 18].
- 3 Disassemble the remaining items and clean them with a lint-free cloth or tissue, see [Disassembling for cleaning ▶ Page 20].
- 4 Reassemble the balance.

**See also**

- 📖 Assembling the pump ▶ Page 18

**5.3.4 Putting into operation after cleaning**

- 1 Reassemble the balance.
- 2 Check if the terminal is connected to the balance.
- 3 Reconnect the balance to the AC/DC adapter.
- 4 Check the level status, level the balance if necessary.
- 5 Respect the warm-up time specified in the "Technical Data".
- 6 Perform an internal adjustment.
- 7 Perform a routine test according to the internal regulations of your company. METTLER TOLEDO recommends to perform a sensitivity test after cleaning the balance.
- 8 Press **→O←** to zero the balance.
  - ➔ The balance is ready to be used.

**See also**

- 📖 Performing an internal adjustment ▶ Page 14

**6 Technical Data****6.1 General data****Power supply**

AC/DC adapter (model no. FSP060-DHAN3):

Input: 100 – 240 V AC ± 10%, 50 – 60 Hz, 1.8 A

Output: 12 V DC, 5 A, LPS, SELV

AC/DC adapter (model no. FSP060-DIBAN2):

Cable for AC/DC adapter:

Balance power consumption:

Polarity:

Input: 100 – 240 V AC  $\pm$  10%, 50 – 60 Hz, 1.5 A

Output: 12 V DC, 5 A, LPS, SELV

3-core, with country-specific plug

12 V DC  $\pm$  10%, 2.25 A

### Protection and standards

Overvoltage category:

II

Degree of pollution:

2

Standards for safety and EMC:

See Declaration of Conformity

Range of application:

Use only indoors in dry locations

### Environmental conditions

The limit values apply when the balance is used under the following environmental conditions:

Height above mean sea level:

Up to 5000 m

Ambient temperature:

+10 – +30 °C

Temperature change, max.:

5 °C/h

Relative air humidity:

30 – 70%, non-condensing

Acclimatization time:

At least **8 hours** after placing the instrument in the same location where it will be put into operation.

Warm-up time:

At least **120 minutes** after connecting the balance to the power supply. When switched on from standby, the instrument is ready for operation immediately.

The balance can be used under the following environmental conditions. However, the weighing performances of the balance may be outside the limit values:

Ambient temperature:

+5 – +40 °C

Relative air humidity:

20% to max. 80% at 31 °C, decreasing linearly to 50% at 40 °C, non-condensing

The balance can be disconnected and stored in its packaging under the following conditions:

Ambient temperature:

-25 – +70 °C

Relative air humidity:

10 – 90%, non-condensing

## 7 Disposal

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties, the content of this regulation must also be related.

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© Mettler-Toledo GmbH 05/2021